TAP SIEVE SHAKER AS 200 TAP

The RETSCH AS 200 tap is suitable for dry sieving with test sieves of 200 mm or 8" diameter. The combination of horizontal, circular sieving motions with vertical taps reproduces the principle of hand sieving. The uniform mechanical action ensures reliable and reproducible measurement results.

This special type of sieving motion used by the AS 200 tap is specified in various standards for particle size analysis of materials such as activated carbon, diamonds, spices, metal powders, abrasives or cement. Operating the AS 200 tap is exceptionally easy and safe. The integrated clamping device allows for sieve stacks with up to 7 or 13 fractions, depending on the height of the sieve frame. The sieving time is set from 1 to 99 minutes via a digital display. The number of rotations and taps is fixed; the tapping motion can be deactivated, if required. A safety switch and an anti-trap protection provide maximum safety. Thanks to an integrated interface, the AS 200 tap can be controlled with the evaluation software EasySieve®.

TAP SIEVE SHAKER AS 200 TAP

AKSESUARLAR VE OPSİYONLAR

RETSCH sieve shakers can be equipped with a variety of accessories to meet a wide range of application requirements.

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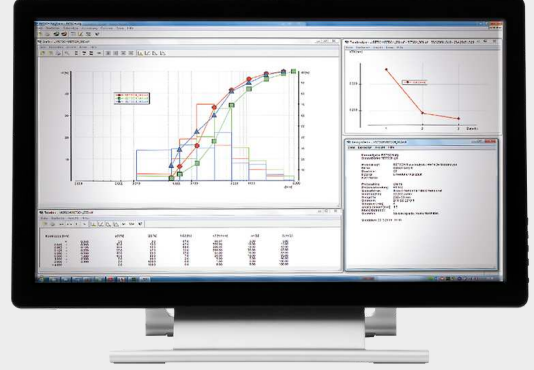
RETSCH TEST SIEVES AND ACCESSORIES - ENGINEERED FOR SUPERIOR PERFORMANCE

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TAP SIEVE SHAKER AS 200 TAP

TIPIK ÖRNEK MALZEMELER

RETSCH's Analytical Sieve Shaker AS 200 tap is ideally suited for separation, fractioning and particle size determination of abrasives, charcoal, cement, coal, diamonds and spices.

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